

### **Importing the Pancancer Pathways Analysis Module**

1. Save the PanCancerPathway1.0.48.zip provided by NanoString to your computer. **Do not** extract the compressed files. Launch nSolver2.5 application.
2. From the “Analysis” menu at the top of the home screen, select “Advanced Analysis Manager”. In the Advanced Analysis Manager window, click “Import New Advanced Analysis” and browse to the directory where the PancancerPathways.zip is saved and click “OK”.
3. Wait for import to complete. The name of the imported Advanced analysis module populates in the Advanced Analysis Manager window. Click “OK” to exit the Advanced Analysis Manager.