

nCounter® Advanced Analysis Installation Note

1. Save the nCounter® Advanced Analysis .ZIP file to your computer. **Do not extract the compressed files.**
2. Launch nSolver™ Analysis Software.
3. From the “Analysis” menu at the top of the home screen, select “Advanced Analysis Manager”.
4. In the Advanced Analysis Manager window, click “Import New Advanced Analysis.”
5. Browse to the directory where the nCounter® Advanced Analysis .ZIP file is saved and click “OK”.
6. Wait for import to complete and the name to populate automatically.
7. Click “OK” to exit the Advanced Analysis Manager.

FOR RESEARCH USE ONLY. Not for use in diagnostic purposes.